



EQUIPMENT REQUIRED:

pH Meter
100 ml beaker
pH 7.00 buffer solution
pH 4.00 buffer solution

PROCEDURE:

1. The pH meter must be calibrated every morning at pH 4.00 and 7.00.
2. To read pH on wine, juice and concentrate, place about 60 ml of sample in a clean 100 ml beaker. Place the pH probe in the beaker and press the measure button. Wait for the meter to display "Ready". Then read the pH meter digital screen and record.

81 M Street
Fresno, CA 93721
Tel: 559-485-2692
Fax: 559-485-4254

640-D Airpark Rd.
Napa, CA 94558
Tel: 707-224-7903
Fax: 707-255-2019

1165 Globe Ave.
Mountainside, NJ 07092
Tel: 908-301-1811
Fax: 908-301-1812

1401 Ware St.
Waupaca, WI 54981
Tel: 715-258-5525
Fax: 715-258-8488