# MLB242 analyser by CDR – Short Guide

Analysis systems based on the MLB242 analyser are:

**CDR FoodLab Junior, CDR WineLab Junior, CDR BeerLab Junior, CDR CiderLab Junior, CDR OxiTester Junior**

**These short instructions do not replace the complete user's manual.**


For the installation of the instrument proceed as follows:

**A.** Place the **MLB242 analyser** on a plane surface in such a position that it is not exposed by mechanical vibrations or extreme temperature fluctuations or excessive light.

**B.** Connect the external power adapter to the port at the rear of the instrument using the appropriate power cord to the mains (110V to 230V).

**C.** After power ON the analyser, the photometer optics will be controlled together with other functional parts.

Appears on the screen:

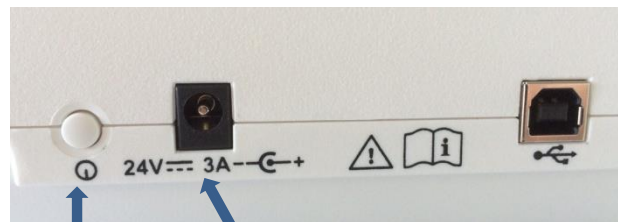ON/OFF button  
To switch OFF press  
for 3 seconds

24V DC electric power port

Test applications are loaded into the working memory for quick access. **1 minute required.**

Followed automatically by next dialogue:

PLEASE WAIT is shown while power up to operating temperature.

**10 minute required.**

After that the **Main dialogue** is shown:

Use "SETTINGS" to configure your system.

Use "PERFORMD TESTS" to get information from Data Base.

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**WARNING!**

Use the device by strictly following the rules and requirements listed below

- Before using the device, all operators should be familiar with the instructions for use that can be downloaded from CDR website.
- Before the first operation, it is recommended that the packaging and the device be visually checked to verify that it has not been damaged during transport. Return the device to the Maintenance department for a functional inspection if damage is reported.
- The device is designed to perform food tests and can be used with cuvettes, code CDR 230012, only.
- The device cannot be used for tests on biological matrices that differ from those described in methods and for which the device has been specifically calibrated during manufacturing.
- The device cannot be used with assay samples that have not been specifically released by CDR.
- The device cannot be used to perform continuous diagnostic tests.
  
- Do not place reflecting objects inside the wells, nor get close to the well opening with your eyes to avoid exposure to light radiations, which are considered, on a prudential basis, as potentially hazardous to the sight.
- The manufacturer recommends to preserve the original box in the event it is necessary to transport the device and install it in another location. In terms of disposal, cardboard is not regarded hazardous for people, animals or materials. Before disposing of the package, see Directive 94/62/EC on packaging and packaging waste.

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CDR S.r.l. Via degli Artigiani, 6